PN: H-020-000168-00

MegaBOLT

生信分析加速器 *Bioinformatics Analysis Accelerator*

快速上手指南
Quick Start Guide

邮箱/E-mail: MGI-service@mgi-tech.com

网址/Website: <https://www.mgi-tech.com/>

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Main Building and Second floor of No.11 Building, Beishan Industrial Zone, Yantian District,
518083 Shenzhen, Guangdong, P.R.China

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---This page is intentionally left blank.---

About this guide

This guide is applicable to MegaBOLT Bioinformatics Analysis Accelerator. The edition is 4.0 and the software version is V2.3.0.0.

This guide and the information contained within are proprietary to MGI Tech Co., Ltd. (hereinafter called MGI), and are intended solely for the contractual use of its customer in connection with the use of the product described herein and for no other purpose. Any person or organization can not entirely or partially reprint, copy, revise, distribute or disclose to others the guide without the prior written consent of MGI. Any unauthorized person should not use this guide.

MGI does not make any promise of this guide, including (but not limited to) any commercial of special purpose and any reasonable implied guarantee. MGI has taken measures to guarantee the correctness of this guide. However, MGI is not responsible for any mistakes or missing parts in the guide, and reserves the right to revise the guide and the software, so as to improve the reliability, performance or design.

Figures in this guide are all illustrations. The contents might be slightly different from the software, please refer to the software purchased.

©2019-2020 MGI Tech Co., Ltd. All rights reserved.

Release date: July 1st, 2021

Manufacturer information

Manufacturer	MGI Tech Co., Ltd.
Manufacturer address	Main Building and Second floor of No.11 Building, Beishan Industrial Zone, Yantian District,518083 Shenzhen, Guangdong, P.R.China
Technical support provider	MGI Tech Co., Ltd.
Technical Support telephone	4000-966-988
Technical support E-mail	MGI-service@mgi-tech.com

Revision history

	Date	Version
Revision	July 1st, 2021	4.0
	December 30, 2020	3.0
	July 17,2020	2.0
Initial release	November 30, 2019	1.0

---This page is intentionally left blank.---

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Contents

1	Quick start	9
	Assembling the device	9
	Logging into the system	9
	Starting the terminal	10
	Viewing Help.....	11
	Submitting a task	11
2	Working with ZLIMS	13
3	Documentation	14
	Brief introduction	14
	Accessing the documents	15
4	Change The Password of Storage Device.....	14

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The image shows a 'Change Passphrase' dialog box with a title bar containing a close button (X). The dialog contains three text input fields: 'Current Passphrase', 'New Passphrase', and 'Confirm Passphrase'. Below the 'Confirm Passphrase' field is a smaller input field with a 'Weak' label to its right. There is an unchecked checkbox labeled 'Show Passphrases'. At the bottom are two buttons: 'Cancel' and 'Change'.

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Assembling the device

Perform the following steps:

1. Connect the server, monitor, keyboard, and mouse.
2. Insert the dongle to a USB port.
3. Connect to the power source and power on the device.

Figure 1 Dongle

Logging into the system

After you power on the device, the user verification interface is displayed, as shown in the figure below:

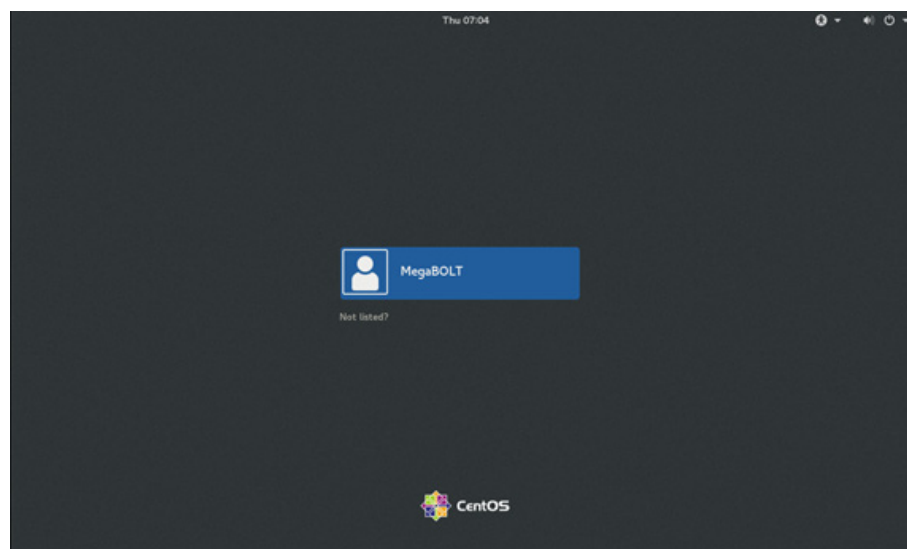
Figure 2 User verification interface

Select the account “MegaBOLT” and input the password “megabolt” to log into the system.

Starting the terminal

To start the terminal, click **Applications>Favorites>Terminal**.

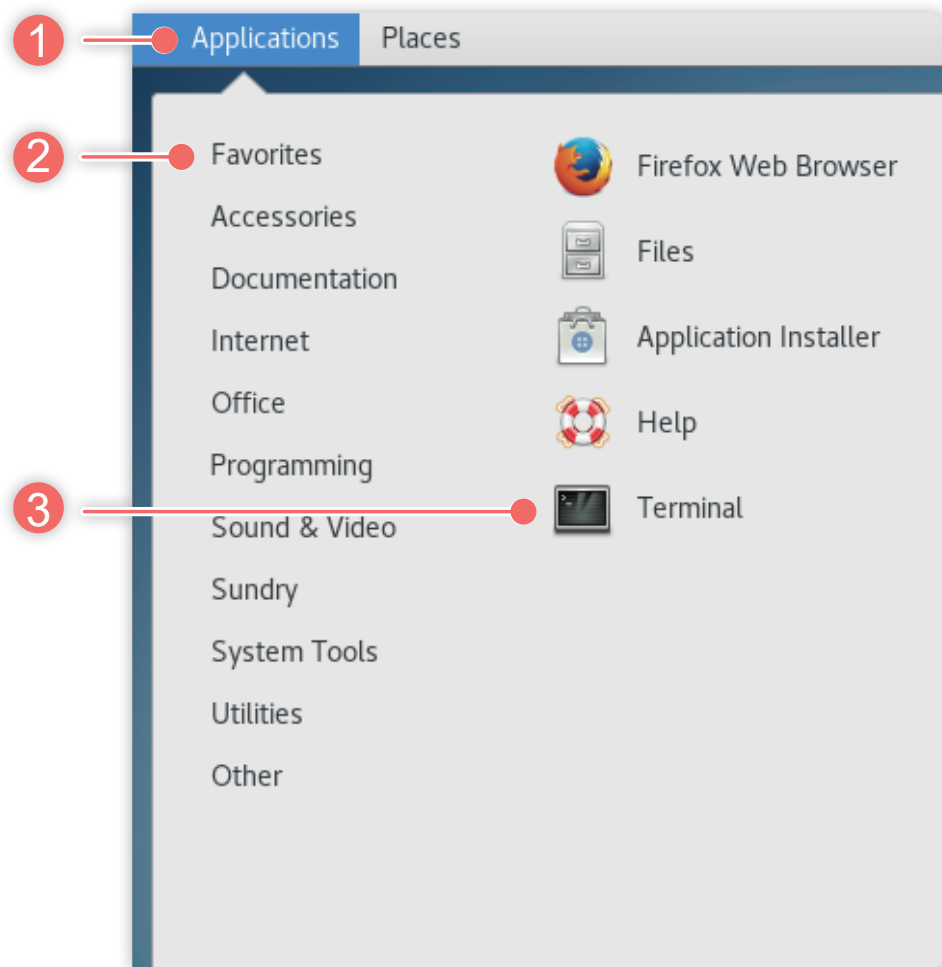
Figure 3 Starting the terminal

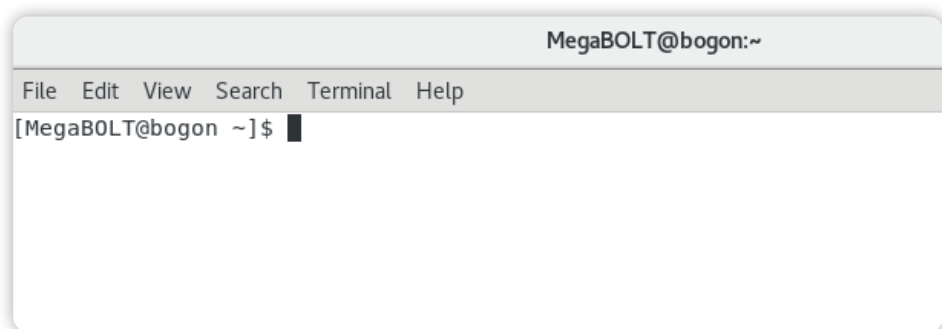
Figure 4 Terminal

Viewing Help

Input the following command on the terminal:

MegaBOLT

Press **Enter** to execute the command and MegaBOLT Help is displayed:

```

MegaBOLT 15:25:41 ~
$ MegaBOLT
Parse argument exception - MegaBOLT help.
MegaBOLT Client Version: v2.0.14.191022_beta-14-gcab686d
Usage: MegaBOLT [Options]
Options:
[common]
--type                <range>  type for pipeline (default: basic)
                                range: <alignment|alignmentsortmarkdup|alignmentsortmarkdupbqsr|
                                bamstats|basic|bqsr|bqsrindex|builddict|buildfai|full|genotype|pegvcfs|
                                haplotypecaller|mutect2|qc|somatic|sortmarkdup|vcfstats>
--help, -h            show help of MegaBOLT
--version, -v         show version of MegaBOLT
--list                string     sample list for SE/PE reads, FORMAT: SampleName Read1_1,Read1_2
                                Read2_1,Read2_2 [adapter1 adapter2] [RGID1,RGID2 RGS1,RGS2
                                RGLB1,RGLB2 RGPL1,RGPL2], no Read2 for SE data
--ref                 string     reference (default: $MEGABOLT_REFERENCE/hg19.fa)
--vcf                 string     dbsnp (default: $MEGABOLT_REFERENCE/dbsnp_151.vcf.gz)
--outputdir, -outdir string     output directory (default: .)
--outputprefix, -prefix string    output file prefix, should not include '/', only used when
                                '--list' is not specified (default: output)
--runtype              <range>  run WGS or WES mode (default: WGS)
                                range: <WES|WGS>
--bed                  <range>  bed file, only used for WES mode (default: BV4)
                                range: <ACV6|ACV6-38|AV2|AV2-38|AV5|AV5-38|AV6|AV6-38|AV7|
                                AV7-38|AIJI|AIJI-38|BV4|BV4-38|BV5|BV5-38|IDT|IDT-38|NME|
                                NME-38|NV3|NV3-38|TV1.2|TV1.2-38|user_defined_path>
                                ACV6      Agilent_V6COSMIC
                                AV2       Agilent_Exome_V2
                                AV5       Agilent_Exome_V5
                                AV6       Agilent_Exome_V6
                                AV7       Agilent_Exome_V7
                                AIJI      AIJI_Exome
                                BV4       BGI_Exome_V4_kit
                                BV5       BGI_Exome_V5_kit
                                IDT       xgen_target
                                NME       Nimblegen_MedExome_V2
                                NV3       Roche_SeqCapEZ_Exome_v3.0
                                TV1.2     Illumina.truseq.v1.2
notice:
1. "*"-"38" is used for hg38, e.g. "BV4-38";
2. user defined bed file path is supported;
do not run markdup in sortmarkdup (default: 0)

```

Figure 5 MegaBOLT Help

For details about the usage of MegaBOLT, refer to *MegaBOLT_User_Manual*.

Submitting a task

MegaBOLT adopts a Server-Client architecture. Analysis tasks are submitted to the MegaBOLT server for unified execution.

1. Execute the following command to run a DEMO analysis task to MegaBOLT:

```
MegaBOLT --list /mnt/ssd/MegaBOLT/test/sample.list
```

The list file contains the sample information to be analyzed, a sample file is shown as below:

```
Test /mnt/ssd/MegaBOLT/test/read1.fq.gz /mnt/ssd/
MegaBOLT/test/read2.fq.gz
```

For detailed format of the *Sample.list* file, refer to *MegaBOLT_User_Manual*.

2. After the analysis task finishes, a folder named after the sample is generated under the current directory to store analysis results.
3. Execute the following command to open the analysis result folder, taking the sample named “Test” for example:

```
cd Test
```

4. Execute the following command to view the result files of the analysis workflow:

```
ll
```

```
Oct 11 20:15 Test.list
Oct 11 20:51 Test.log
Oct 11 20:42 Test.mm2.sortdup.bqsr.bam
Oct 11 20:42 Test.mm2.sortdup.bqsr.bam.bai
Oct 11 20:42 Test.mm2.sortdup.bqsr.bam.grp
Oct 11 20:51 Test.mm2.sortdup.bqsr.HaplotypeCaller.vcf.gz
Oct 11 20:51 Test.mm2.sortdup.bqsr.HaplotypeCaller.vcf.gz.out
Oct 11 20:51 Test.mm2.sortdup.bqsr.HaplotypeCaller.vcf.gz.tbi
Oct 11 20:51 Test.out
```

Figure 6 Result files

If you have any doubts about the analysis workflow, refer to *MegaBOLT_User_Manual* or contact the technical support.

2 Working with ZLIMS

You can log into ZLIMS (Zebra Laboratory Information Management System) to assign a MegaBOLT analysis task and check the analysis report. ZLIMS adopts an easy-to-use web design. You can submit tasks and check the results on the website easily even if you have no bioinformatics background. For details, refer to *MGI ZLIMS (Lite) User Manual__Standard V2.0.5*.

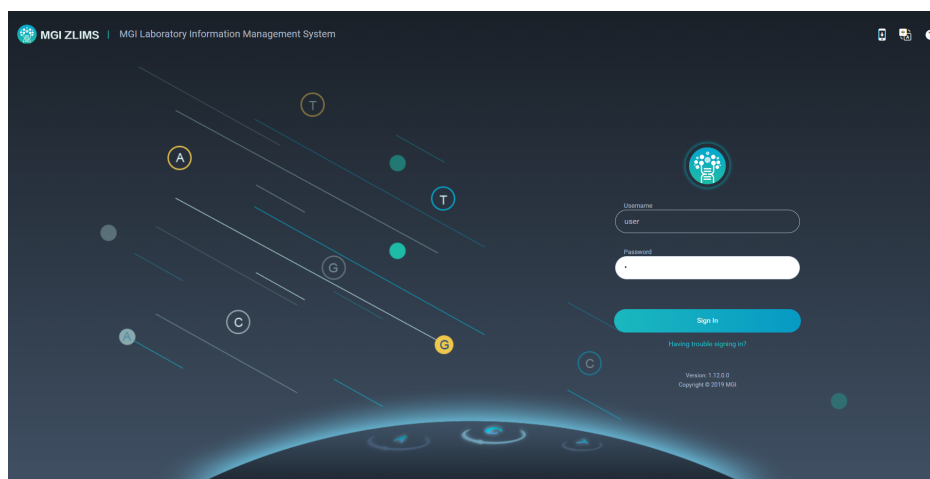
Figure 7 ZLIMS login page

3

Documentation

Brief introduction

The following documentation is provided with MegaBOLT:

Table 2 MegaBOLT documentation

Name	Description
MegaBOLT_Quick_Start_Guide	It is intended to guide the user through task analysis workflow and result confirmation, and provide basic information for the other three files.
MegaBOLT_User_Manual	It provides information about MegaBOLT's basic features and its functions with samples.
MegaBOLT_User_Manual_ZLIMS	ZLIMS is a laboratory management system of MGI. You can log into ZLIMS to assign a MegaBOLT analysis task and check the analysis results.
MegaBOLT_User_Manual_Advanced	It provides information about the storage structure, storage mounting, source code path for some software tools, and FAQ. It further guides the user to develop workflow with MegaBOLT and its submodules.

Accessing the documents

A digital version of this guide is available on the desktop. You can find the other documents by clicking **Home>Documents** on the desktop, as shown in the figure below:

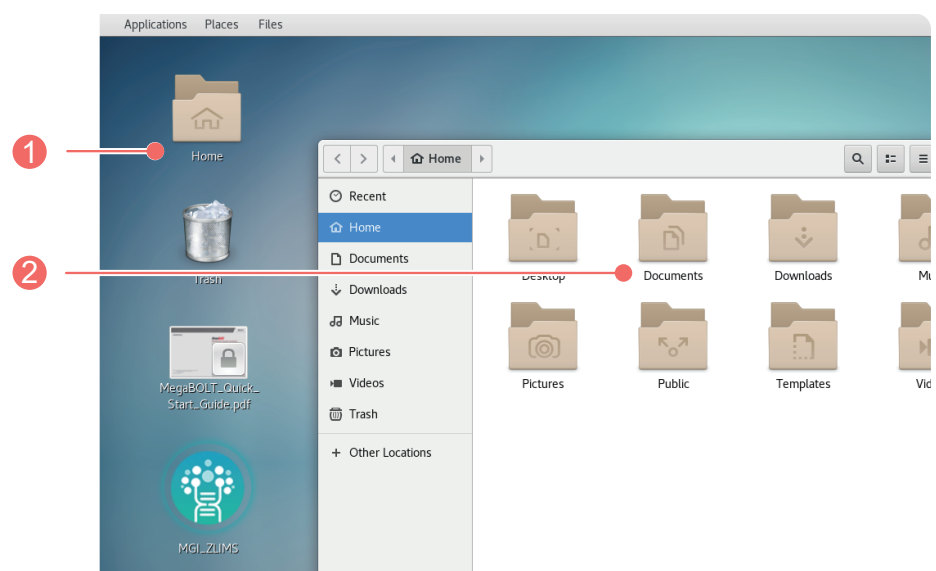
Figure 8 Document directory

You can also obtain the documents from the following directory:

/home/MegaBOLT/Documents

4

Change The Password of Storage Device

For regions where GDPR (General Data Protection Regulation) is applied, please change the initial passphrase on your first entrance to MegaBOLT system, and perform the following processes. For more details about data storage device encryption and decryption, please refer to Chapter 10 of MegaBOLT_User_Manual_Advanced.

1. Open the Linux Disks tool by the following steps:
Desktop → click Applications → click Utilities → click Disks
2. Choose the encrypted disk data, click the setting button, and then click Change Passphrase on the prompted menu.

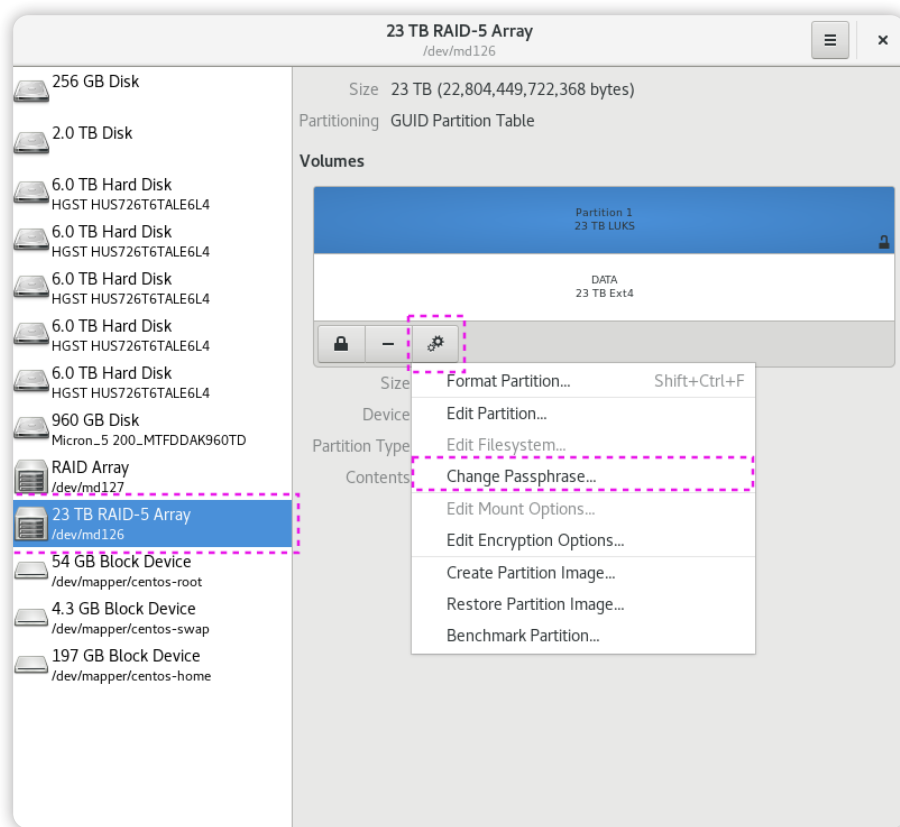
图 9 Change PWD 1

3. Enter your new password and then confirm it, click Change when all setup is ready.

Change Passphrase

Current Passphrase

New Passphrase

Confirm Passphrase

Weak

☐

Show Passphrases

Cancel

Change

图 10 Change PWD 1

Warning

Please keep the new password properly and share it among administrators. Once the password is lost, all data in the encrypted storage device cannot be recovered. You have to reformat the device before you use it again.